# CAMCOLMS4

**MOBILE PHONE MICROSCOPE  
 MICROSCOOP VOOR SMARTPHONE  
 MICROSCOPE POUR SMARTPHONE  
 MICROSCOPIO PARA SMARTPHONE  
 MIKROSKOP FÜR SMARTPHONE  
 MIKROSKOP DO TELEFONÓW KOMÓRKOWYCH  
 MICROSCÓPIO PARA TELEMÓVEL**

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# USER MANUAL

## 1. Introduction

To all residents of the European Union

### Important environmental information about this product

This symbol on the device or the package indicates that disposal of the device after its lifecycle could harm the environment. Do not dispose of the unit (or batteries) as unsorted municipal waste; it should be taken to a specialized company for recycling. This device should be returned to your distributor or to a local recycling service.

Respect the local environmental rules.

### If in doubt, contact your local waste disposal authorities.

Thank you for choosing Velleman! Please read the manual thoroughly before bringing this device into service. If the device was damaged in transit, do not install or use it and contact your dealer.

## 2. Safety Instructions

- Keep this device away from children and unauthorized users.
- This device can be used by children aged from 8 years and above, and persons with reduced physical, sensory or mental capabilities or lack of experience and knowledge if they have been given supervision or instruction concerning the use of the device in a safe way and understand the hazards involved. Children shall not play with the device. Cleaning and user maintenance shall not be made by children without supervision.
- Children shall be supervised to ensure that they do not play with the appliance.
- This device contains small parts. Choking hazard!
- The white LED, which illuminates the microscope target field, is very bright. Do not stare directly into this LED as it may damage your eyes.

## 3. General Guidelines

Refer to the Velleman® Service and Quality Warranty on the last pages of this manual.

- Keep this device away from dust and extreme temperatures.

- Protect this device from shocks and abuse. Avoid brute force when operating the device.

- Familiarise yourself with the functions of the device before actually using it.
- All modifications of the device are forbidden for safety reasons. Damage caused by user modifications to the device is not covered by the warranty.
- Only use the device for its intended purpose. Using the device in an unauthorised way will void the warranty.
- Damage caused by disregard of certain guidelines in this manual is not covered by the warranty and the dealer will not accept responsibility for any ensuing defects or problems.
- Do not switch the device on immediately after it has been exposed to changes in temperature. Protect the device against damage by leaving it switched off until it has reached room temperature.
- Keep this manual for future reference.

## 4. Overview

Refer to the illustrations on page 2 of this manual.

<b>1</b>	body
<b>2</b>	zoom dial
<b>3</b>	turret
<b>4</b>	objectives
<b>5</b>	slide clips
<b>6</b>	base LED
<b>7</b>	specimen stage

<b>8</b>	base
<b>9</b>	arm
<b>10</b>	focusing knob
<b>11</b>	reflective LED
<b>12</b>	storage drawer
<b>13</b>	LED on-off switch

## 5. Battery

### WARNING

Do not puncture batteries or throw them in fire as they may explode. Do not attempt to recharge non-rechargeable batteries (alkaline). Dispose of batteries in accordance with local regulations. Keep batteries away from children.

1. Open the battery compartment.
2. Install three fresh AA batteries following the polarity markings inside the compartment.
3. Close the battery compartment.
4. Remove the batteries from the device when not in use.

## 6. Operation

1. Place a prepared specimen slide on the specimen stage and slide it into the clips, such that the clips hold the slide securely.
2. Switch on the LED – the base LED or reflective LED – to illuminate the specimen.
3. Turn the objective turret – the turret can be rotated both ways – to the desired magnification. Rotate the objective turret such that the objective is directly on top of the specimen. It is recommended to start with the least powerful magnification rate and select a more powerful objective if needed.
4. View through the eyepiece and rotate the focusing knob clockwise or anticlockwise until the image is sharp. Make sure the objective is clear from obstacles and does not hit the specimen while lowering the turret.
5. Zoom in on the image with the zoom dial.

### REMARK

The arm can be swivelled for more ease-of-use.  
The body can be removed from the arm for transportation.  
Viewing through the eyepiece requires sufficient light.

## 7. Using the Microscope with your Smartphone

You can easily take pictures of your specimen using your smartphone. To do so, slide the smartphone platform onto the body and place the smartphone onto the platform in such way, that the camera sits just above the eyepiece.

Now you can take pictures, record and even zoom through your smartphone.

## 8. Cleaning and Maintenance

Occasionally wipe with a damp cloth to keep it looking new. Do not use harsh chemicals, cleaning solvents or strong detergents.

## 9. Technical Specifications

magnification rate .....	50, 100, 200, 400x
battery .....	3 x LR6/AA battery (not incl.)
weight.....	312 g
dimensions .....	96 x 135 x 225 mm

**Use this device with original accessories only. Velleman nv cannot be held responsible in the event of damage or injury resulting from (incorrect) use of this device. For more info concerning this product and the latest version of this manual, please visit our website [www.velleman.eu](http://www.velleman.eu). The information in this manual is subject to change without prior notice.**

### © COPYRIGHT NOTICE

**The copyright to this manual is owned by Velleman nv. All worldwide rights reserved.** No part of this manual may be copied, reproduced, translated or reduced to any electronic medium or otherwise without the prior written consent of the copyright holder.

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**EN****Velleman® Service and Quality Warranty**

Since its foundation in 1972, Velleman® acquired extensive experience in the electronics world and currently distributes its products in over 85 countries.

All our products fulfil strict quality requirements and legal stipulations in the EU. In order to ensure the quality, our products regularly go through an extra quality check, both by an internal quality department and by specialized external organisations. If, all precautionary measures notwithstanding, problems should occur, please make appeal to our warranty (see guarantee conditions).

**General Warranty Conditions Concerning Consumer Products (for EU):**

- All consumer products are subject to a 24-month warranty on production flaws and defective material as from the original date of purchase.
- Velleman® can decide to replace an article with an equivalent article, or to refund the retail value totally or partially when the complaint is valid and a free repair or replacement of the article is impossible, or if the expenses are out of proportion. You will be delivered a replacing article or a refund at the value of 100% of the purchase price in case of a flaw occurred in the first year after the date of purchase and delivery, or a replacing article at 50% of the purchase price or a refund at the value of 50% of the retail value in case of a flaw occurred in the second year after the date of purchase and delivery.

**• Not covered by warranty:**

- all direct or indirect damage caused after delivery to the article (e.g. by oxidation, shocks, falls, dust, dirt, humidity...), and by the article, as well as its contents (e.g. data loss), compensation for loss of profits;
- consumable goods, parts or accessories that are subject to an aging process during normal use, such as batteries (rechargeable, non-rechargeable, built-in or replaceable), lamps, rubber parts, drive belts... (unlimited list);
- flaws resulting from fire, water damage, lightning, accident, natural disaster, etc....;
- flaws caused deliberately, negligently or resulting from improper handling, negligent maintenance, abusive use or use contrary to the manufacturer's instructions;
- damage caused by a commercial, professional or collective use of the article (the warranty validity will be reduced to six (6) months when the article is used professionally);
- damage resulting from an inappropriate packing and shipping of the article;
- all damage caused by modification, repair or alteration performed by a third party without written permission by Velleman®.
- Articles to be repaired must be delivered to your Velleman® dealer, solidly packed (preferably in the original packaging), and be completed with the original receipt of purchase and a clear flaw description.
- Hint: In order to save on cost and time, please reread the manual and check if the flaw is caused by obvious causes prior to presenting the article for repair. Note that returning a non-defective article can also involve handling costs.
- Repairs occurring after warranty expiration are subject to shipping costs.
- The above conditions are without prejudice to all commercial warranties.

**The above enumeration is subject to modification according to the article (see article's manual).**

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